


# BROEN-LAB ESS

## Mounting instruction for products and signs

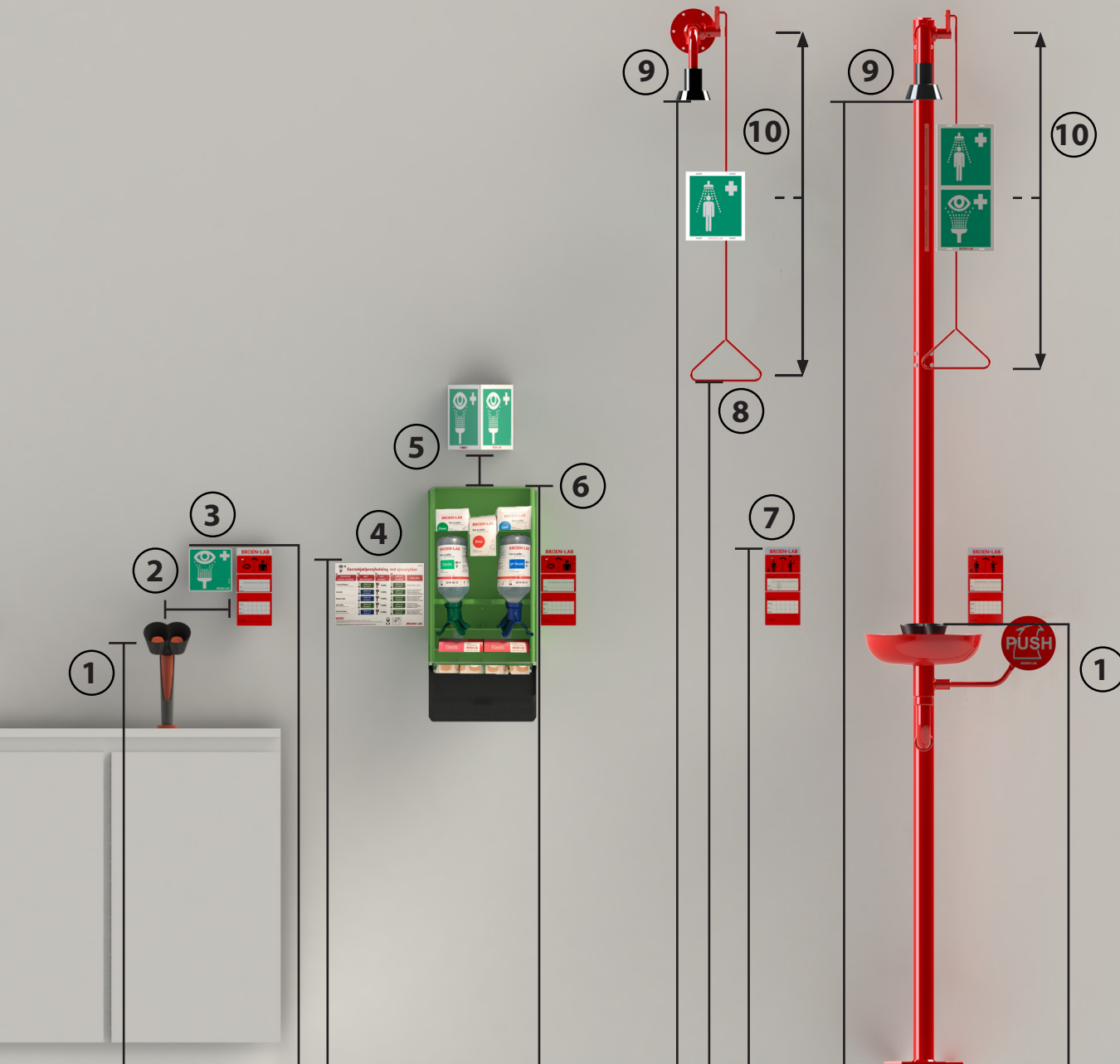

- ① Eye shower: **1000 mm ± 200 mm above the level on which the user stands**
- ② Distance between sign and product: **Max. 250 mm**
- ③ Sign on the wall: **Max. 2000 mm over the floor**
- ④ First aid plan: **Max. 1500 mm over the floor**
- ⑤ Triangular sign: **Min. 250 mm straight above the product**
- ⑥ Eye Wash Station: **Max. 1700 mm over the floor**
- ⑦ Service sign: **Max. 1400 mm over the floor**
- ⑧ Pullrod: **Max. 1750 mm over the floor**
- ⑨ Body shower: **2200 mm ± 100 mm above the level on which the user stands**
- ⑩ Sign placed: **Center of pullrod**

**BROEN-LAB**