

BROEN-LAB service visit

Overview of the contents

All emergency showers are checked according to EN15154 part 1, 2 & 5 as described below.

First service visit

All emergency showers are assessed according to our assessment guide

- Accessibility and tidiness of the area
- Shower location
- Signs and symbols
- Valve function
- Spray image from shower head and eye wash station is tested visually with testing equipment
- Test and check of dust covers
- Test for water flow
- Visual check of water quality
- Cleaning of emergency shower where emergency shower is wiped off for dust and impurities

BROEN-LAB assessment report with photo documentation is filled out and sent to the person in charge of safety at your organization and inspection sticker is filled out and attached to all the inspected showers.

Following service visits

All emergency showers are assessed according to our service report

- Accessibility and tidiness of the area
- Shower location
- Signs and symbols
- Valve function
- Spray image from shower head and eye wash station is tested visually with testing equipment
- Test and check of dust covers
- Test for water flow
- Visual check of water quality
- Cleaning of emergency shower where emergency shower is wiped off for dust and impurities

BROEN-LAB service report with photo documentation is filled out and sent to the person in charge of safety at your organization and inspection sticker is filled out and attached to all the inspected showers.